

Silent Installation Instructions

Instructions on how to create a silent installation file for DinoCapture 2.0 software or device driver of the Dino-Lite Pro series.

Notice: The AM211, AM311, and AM313 series may not be suitable for silent installation.

Steps

1. To start the silent installation process, please download **DNC_SI** from:
www.dino-lite.com/support/DNC2_SI_v1513.exe
2. DNC2_SI.exe can be unpacked by double clicking on it after downloading.
3. Open the Windows MS -DOS command prompt shell window (cmd).
4. In the command prompt, change the directory to where the DNC2_AP is located.
5. Record a response file of the setup.exe in the DNC2_AP folder by typing
`setup.exe /r` in the command prompt.
6. The DinoCapture 2.0 installshield will then start. When finished with the installation, the Silent Response file **setup.iss** will be created in the *C:\Windows* folder.
7. Move the setup.iss file into the DNC2_AP folder.
8. You can now run the DinoCapture 2.0 software installation in silent mode by typing `setup.exe /s` in the command prompt.
9. If plan to create a silent response file for the Dino-Lite Pro series device driver, repeat steps 3 to 8 but change the directory to DN4xxDrv.